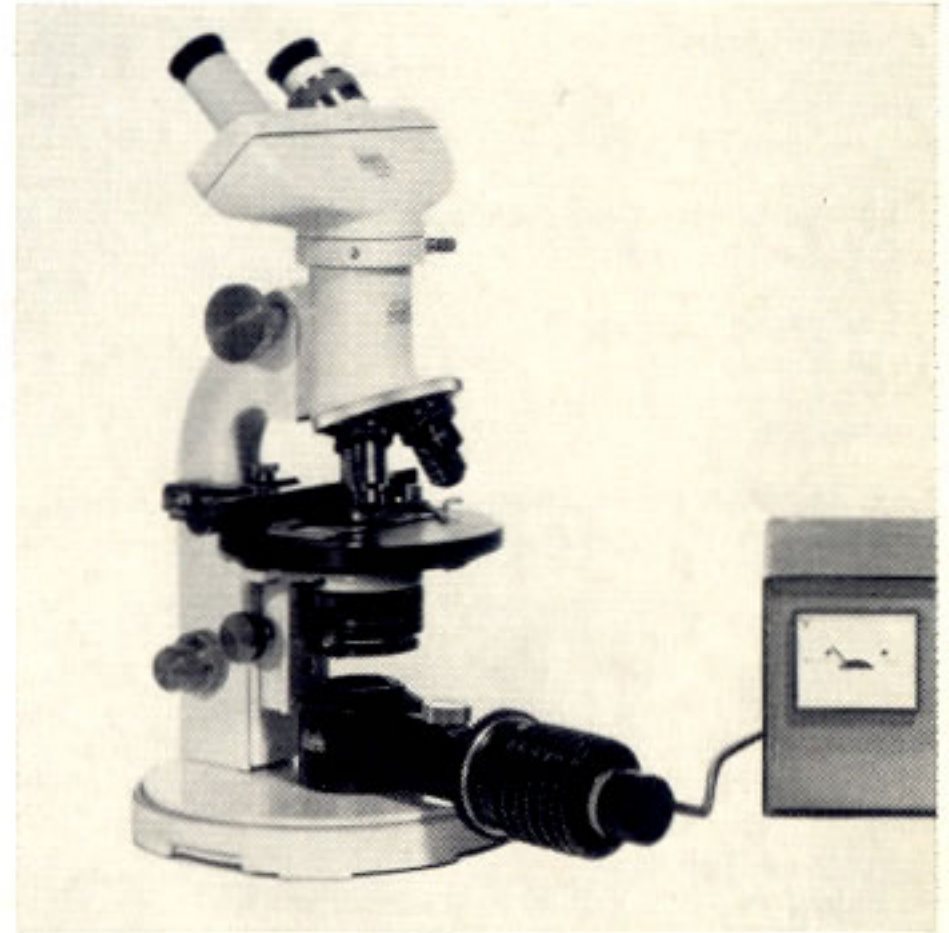
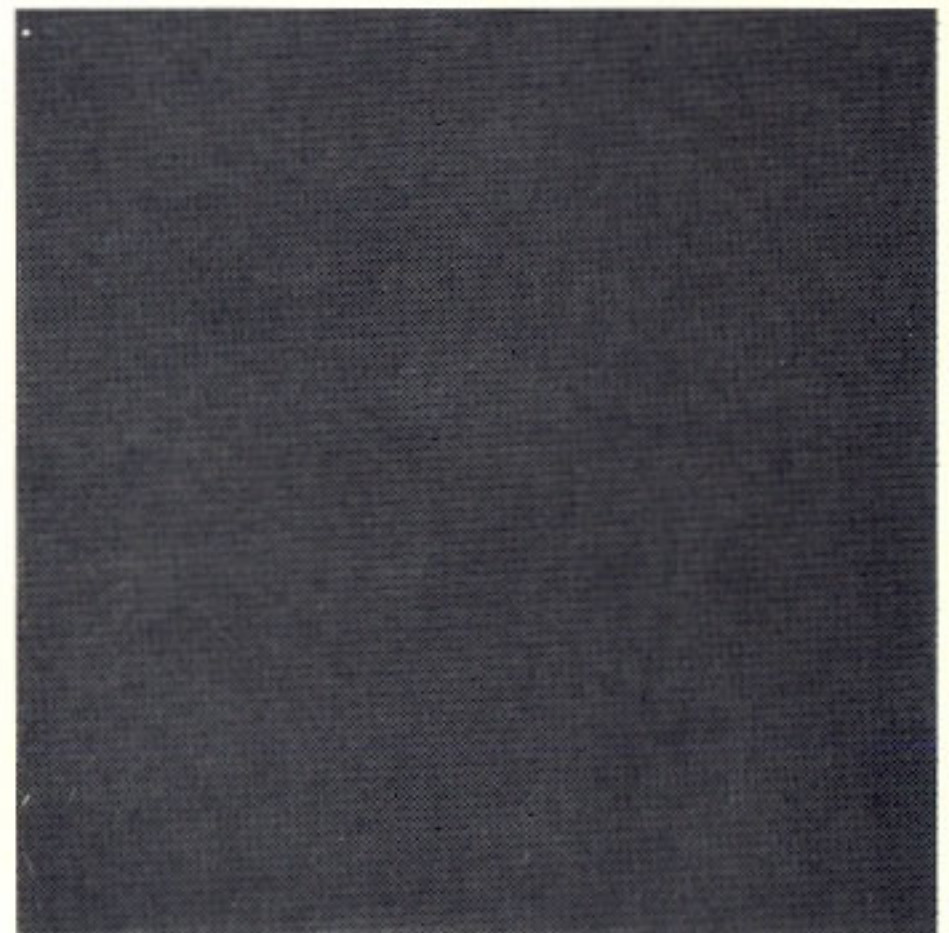
Microscope
Microscope
Mikroskop
Microscopio

WILD M 11

Instructions for use
Mode d'emploi
Gebrauchsanweisung
Instrucciones para el uso

WILD
HEERBRUGG

SWITZERLAND

Contents	Page
Unpacking	2
Assembly	3
Setting up the illumination	7
Mirror	7
Mains (line) lamp M	7
Socket lamp P	8
Transformer	9/10
Attachable Köhler lamp Q	10
Adjusting the binocular tube	12
Care of the microscope	13
Packing	14

The numbers quoted in the text refer to the fold-out illustration on page 16.

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Unpacking

Open catches (a) of metal hood and lift hood off with care. Remove padding, and remove protective covers from tubes. Raise microscope body using coarse focussing control (1).

If a **wooden cabinet** is purchased, first remove the securing screws from the underside.

Important! Free-running of coarse adjustment can be regulated by simultaneously turning the two coarse focussing knobs (1) in opposite directions.

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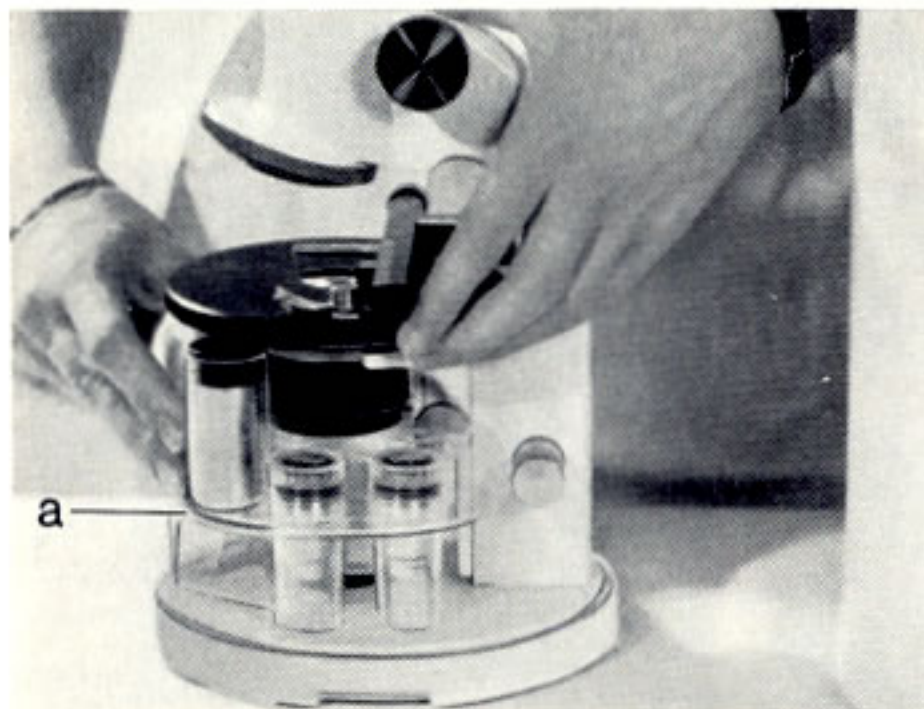
Optics holder / Fine focussing protection

The optics holder (a) has a double function during transport. It stores the objectives and eyepieces and also takes the load off the fine focussing mechanism. Lift stage manually and remove optics holder (a). Screw objectives into nosepiece and fit eyepieces into tubes.

Changing tubes

Monocular tube is screwed on to microscope body, and is rotatable through 360°. It is removed by turning milled ring anticlockwise. To fit binocular tube, an adapter (a) must first be fitted. Screw off monocular tube: screw adapter on to microscope body (b).

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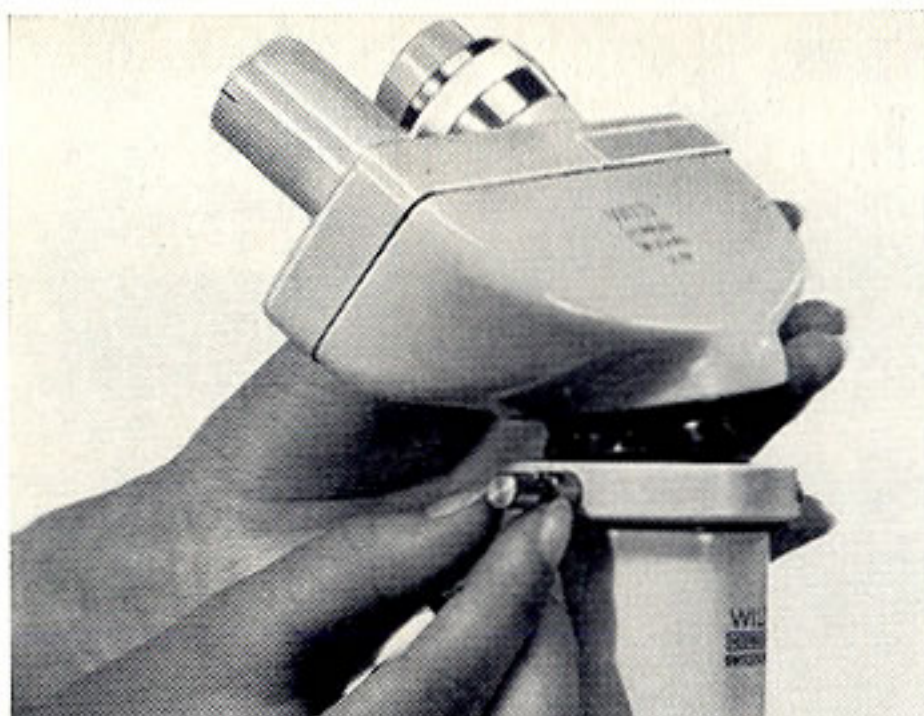
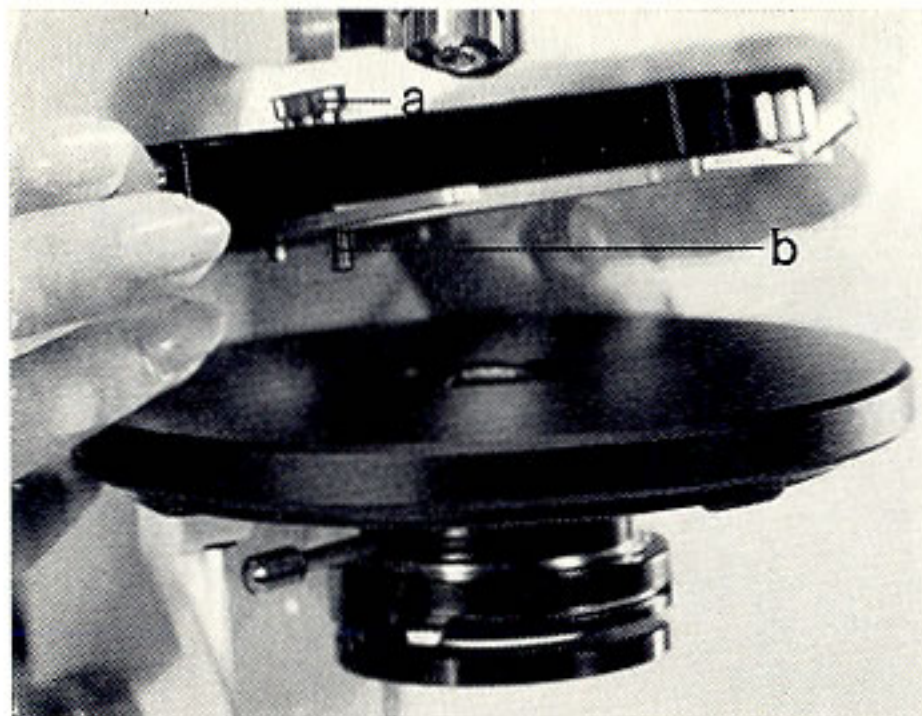
Loosen clamping screw (5) of adapter. Press binocular tube against spring bolt of clamping screw (5) and lower it into position. Turn tube to desired position (rotatable through 360°) and retighten clamping screw.

Fitting mechanical stage

Place mechanical stage on microscope stage so that guide pins (b) of former engage corresponding groove or hole in latter. Mechanical stage is secured by means of screw.

Note: When replacing metal hood, ensure that coordinates of mechanical stage are roughly in the middle of their range of movement.

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
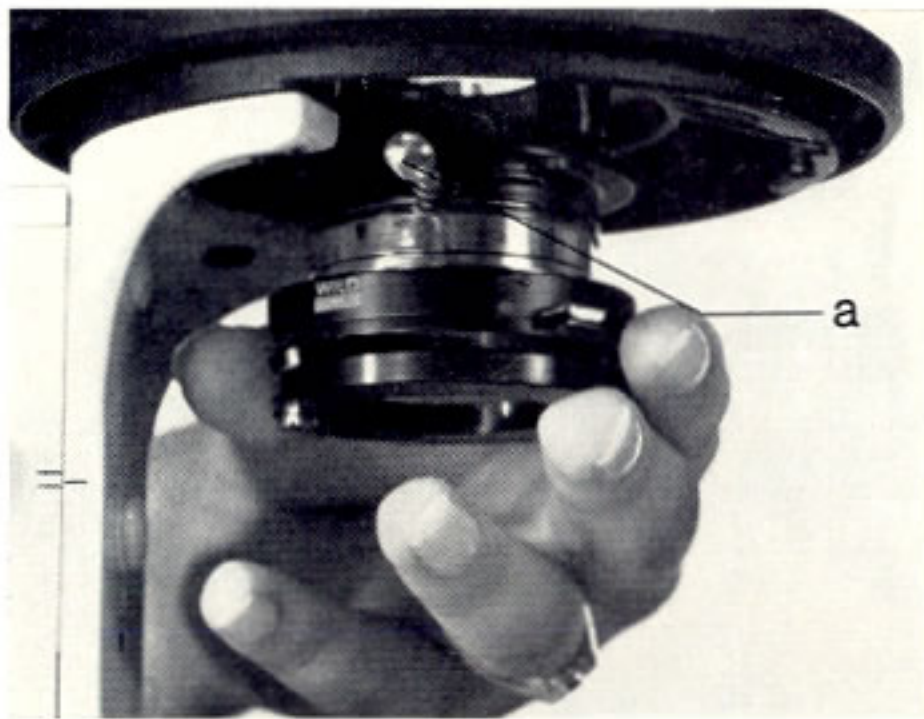
Fitting condenser (Stand A)

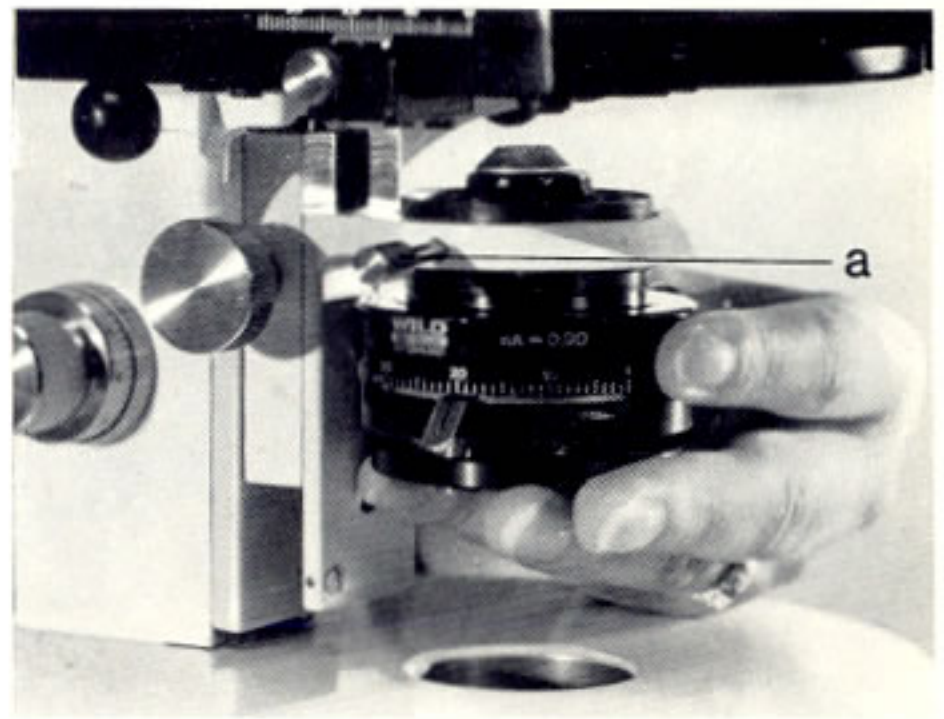Condenser holder of Stand A is immovably fixed beneath microscope stage. Loosen screw (a). Push condenser fully into holder and secure it with clamping screw (a).

Condenser (Stand B)

Lower condenser holder somewhat by means of condenser drive (2). Loosen clamping screw (a). Push condenser fully home and retighten screw.

Note: Front lens of the 0.65/1.30 swing-out condenser must be swung out during fitting: swing-out knob must be diametrically opposite to clamping screw for all condensers. Tighten clamping screw.

Free-running of condenser drive can be regulated by turning its two drive knobs (2) in opposite directions.

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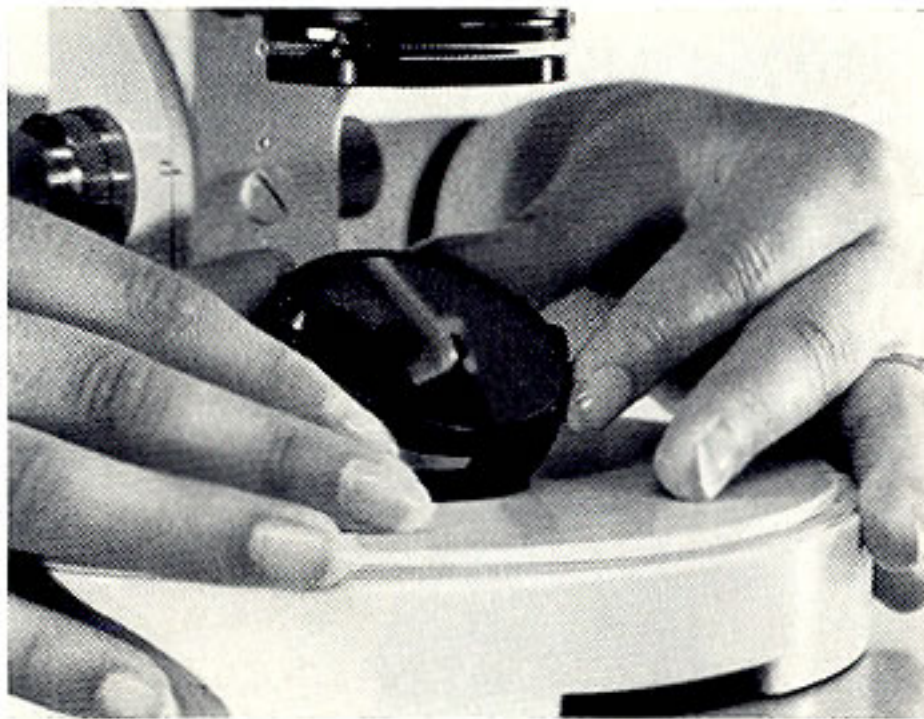
Mirror

Set plane side of mirror to face light. Fully raise condenser by means of condenser drive (2) and open condenser diaphragm (6) completely. Move mirror until field of view through eyepiece appears bright.

Focus specimen sharply with coarse and fine adjustments (1 and 3). Readjust mirror until field of view again appears uniformly illuminated.

If a separate light source is used, bulb filament should be focussed as sharply as possible in middle of plane side of mirror, by adjusting lamp socket or collector. Beyond this, instructions for Köhler illumination are to be followed.

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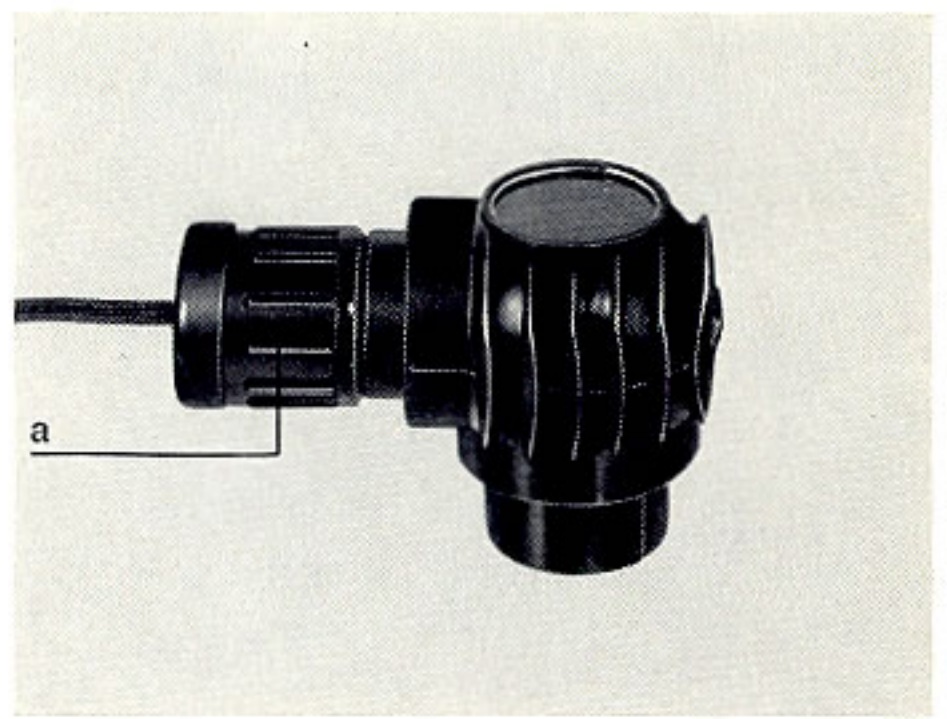
Mains (line) lamp M

Connect lamp to mains (line) supply. Switch on lamp by turning ridged collar (a). Fully raise condenser and fully open condenser diaphragm (6). Focus specimen. Adjust condenser diaphragm according to instructions on page 11.

Changing bulb: Withdraw lamp socket from housing: unscrew and replace bulb. When replacing socket, note guide groove.

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
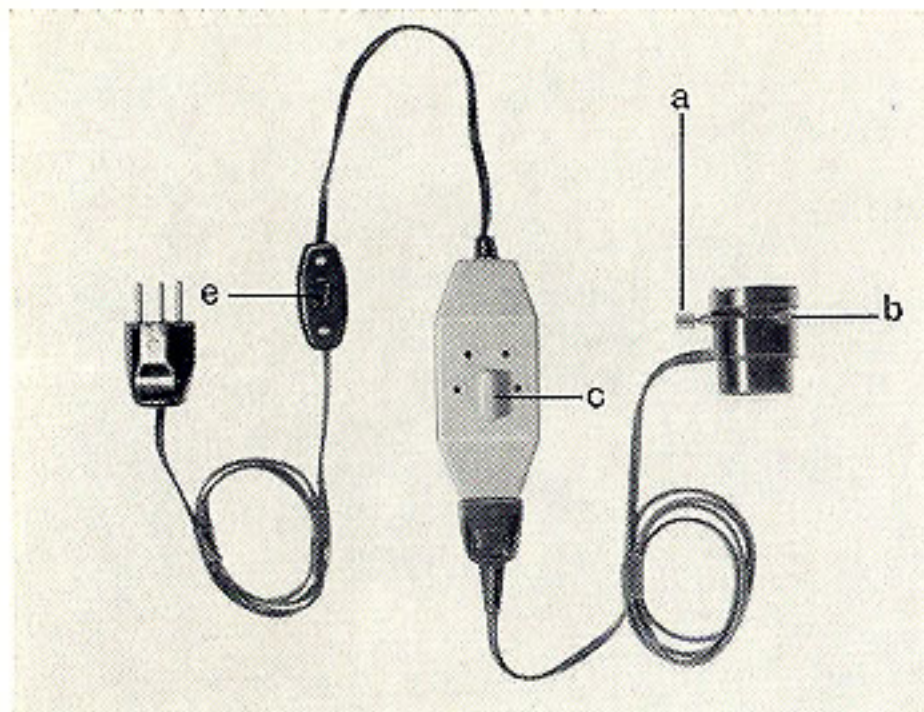
Socket lamp P, 6V/5W

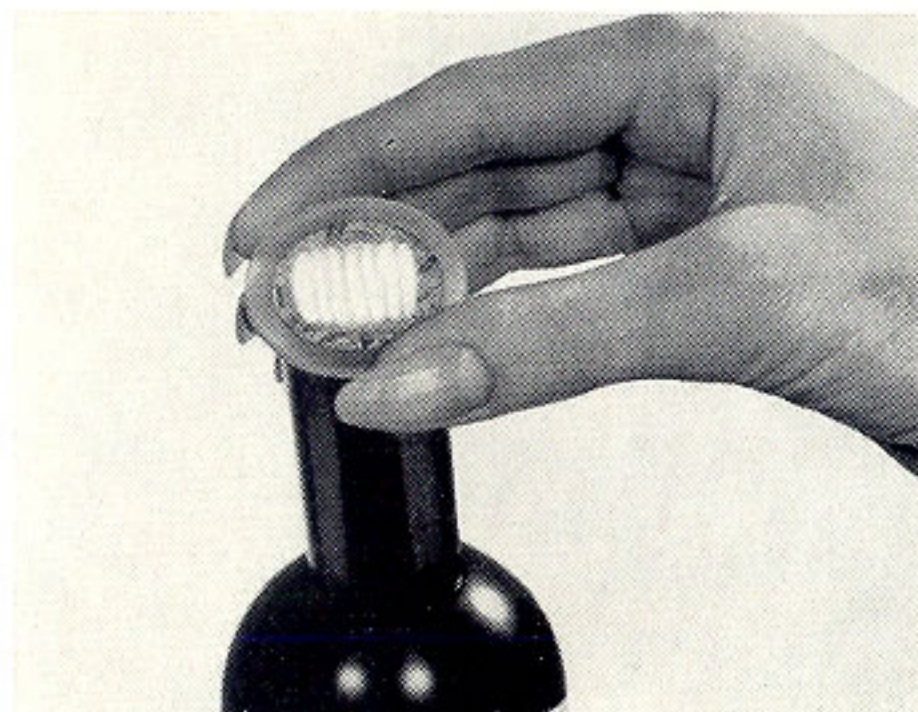Set voltage selector on transformer (c) to mains (line) voltage. Connect lamp to transformer and transformer to mains (line). Switch on (e). Place specimen on stage, fully raise condenser and open condenser diaphragm (6).

Place frosted filter over empty tube. Using a 10 $\times$ objective, focus image of specimen on filter. Lower condenser until lamp filament is imaged on filter. Using the two centring screws (a), centre filament image in tube aperture. Place filter in lamp, insert eyepiece and refocus with fine adjustment. Adjust condenser diaphragm according to instructions on page 11.

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Changing bulb

Loosen both centring screws (a) of socket lamp by about two turns. Press collector mount against counterspring (b) and lift upwards out of lamp housing. Press bulb downwards and release from bayonet socket by means of a slight turn to left. Press new bulb into place and turn it to right. Replace collector mount.

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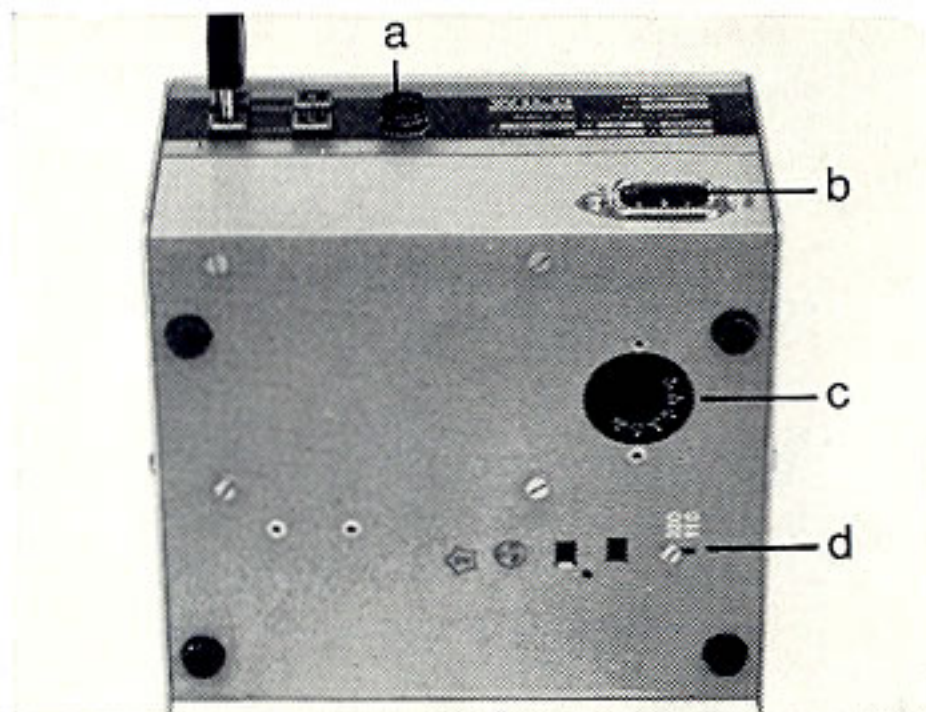
Regulating transformer f. Köhler lamp Q, 6V/20W

Connect cable of lamp to a secondary output of transformer. Set transformer to mains (line) voltage, as follows:

1. Turn voltage selector (c) to correct mains (line) voltage by means of a coin.
2. Loosen screw (d) slightly, slide it to correct voltage setting, and retighten it.

For voltages from 110 to 150V, use position "110". For voltages from 220 to 250V, use position "220". Plug mains (line) cable into socket (b) and connect other end to mains (line). Switch on transformer by turning regulating knob clockwise past click-stop at 2V, and set it to 6V, at which position there is another click-stop. Fuses (a): 6.3A, 6.3 x 32 mm

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
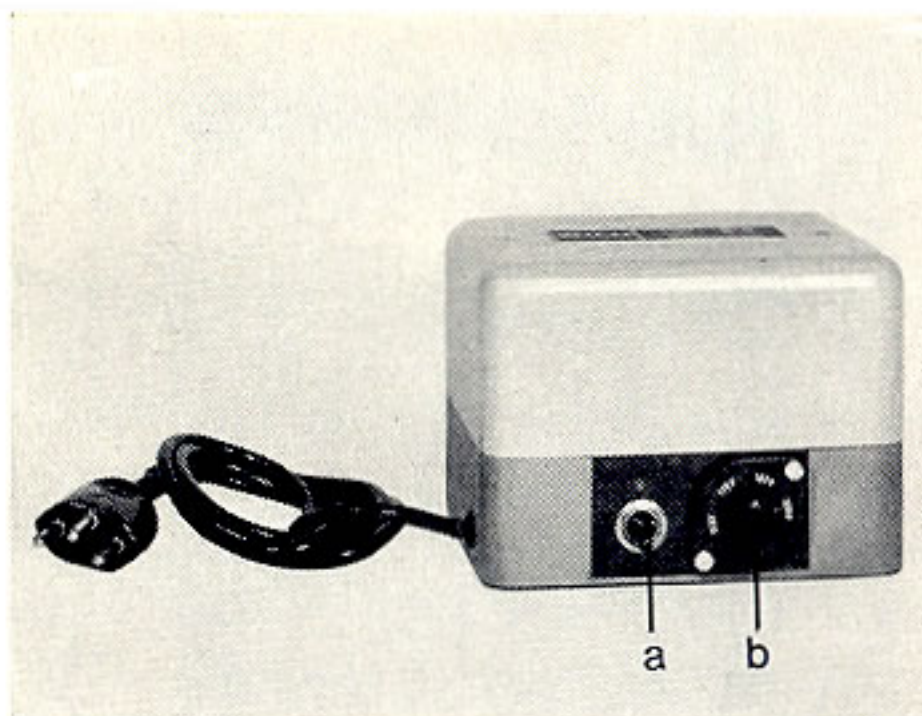
Step transformer for Köhler lamp Q, 6V/20W

Connect cable of lamp to 6V secondary output of transformer. Remove fuse (b) from voltage selector by pressing it in gently and turning it to the left. Set voltage selector to available mains (line) voltage by means of a coin. Press fuse in and secure it by turning it to the right.

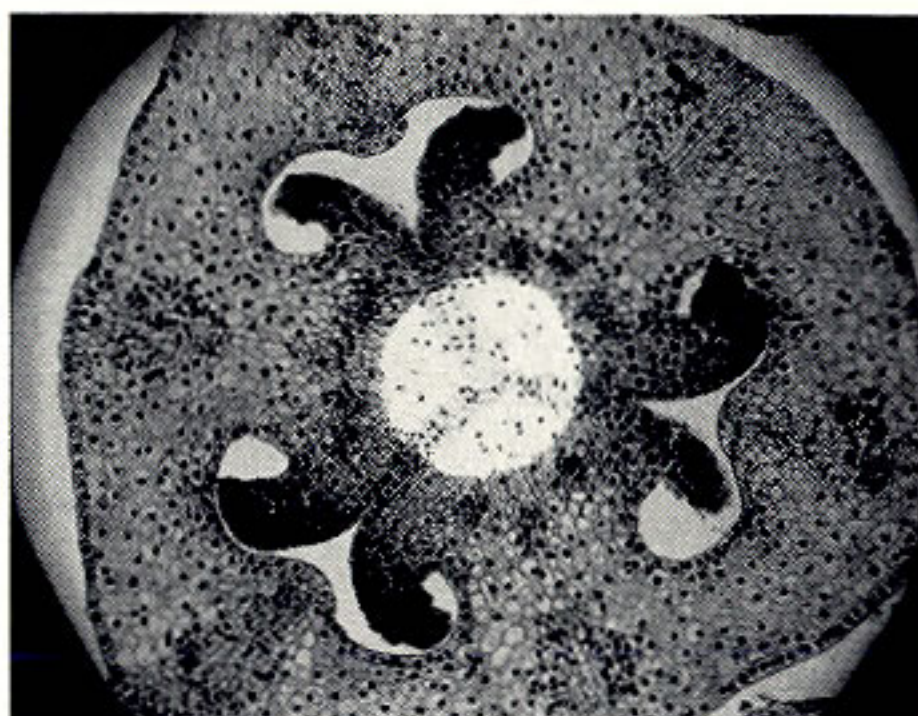Connect mains (line) cable of transformer to mains (line). Switch on transformer by means of switch (a). Fuse: 0.5A, 6.3 x 32 mm

Attachable Köhler lamp Q, 6V/20W

Select 10x objective, raise condenser (with front lens) to highest position and open condenser diaphragm (6). Focus a specimen. Half close field diaphragm (9) and lower condenser until field diaphragm is sharply imaged in specimen plane. Turn and release centring knob (8), centre image of field diaphragm. Reclamp knob (8).

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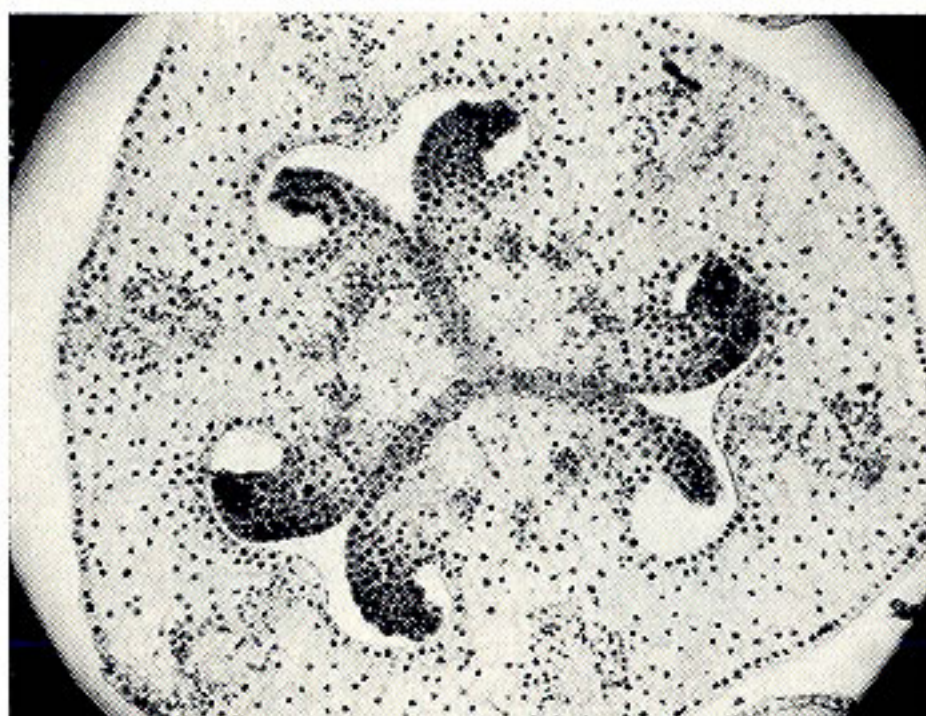
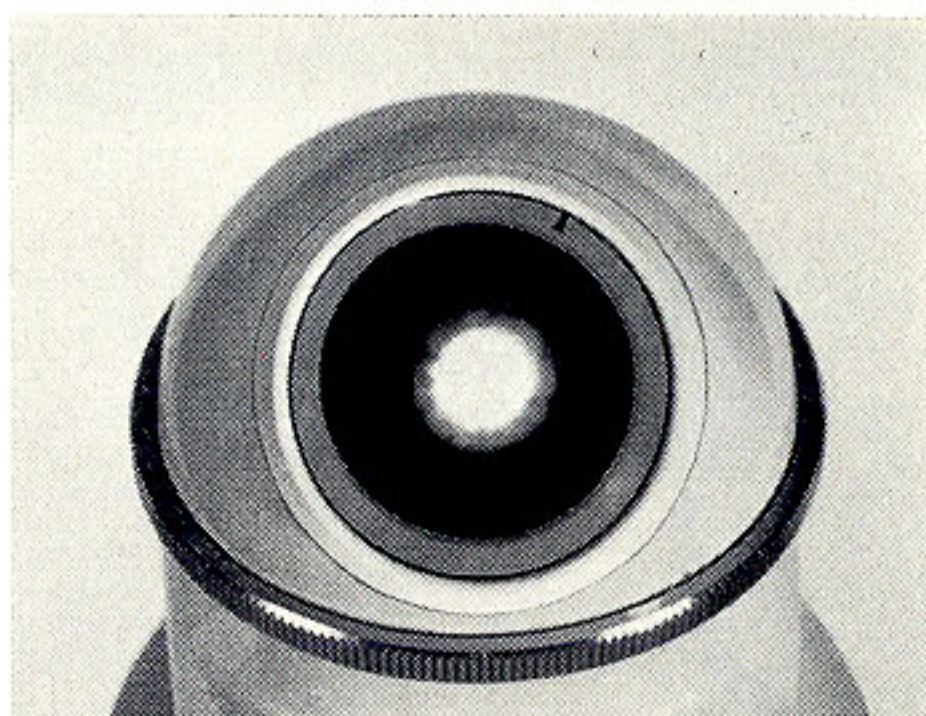
Open field diaphragm (9) until it just disappears from field of view. If it is not possible to illuminate whole field, front lens of the condenser must be removed and diaphragm and condenser adjustments must be repeated.

Adjusting condenser diaphragm

Remove eyepiece and look into empty tube from a distance of about 10 in. (25 cm). Slowly close condenser diaphragm (6) until about 1/3 of area of light spot is covered. Replace eyepiece. Condenser diaphragm and field diaphragm (if fitted) must be readjusted for each change of objective.

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
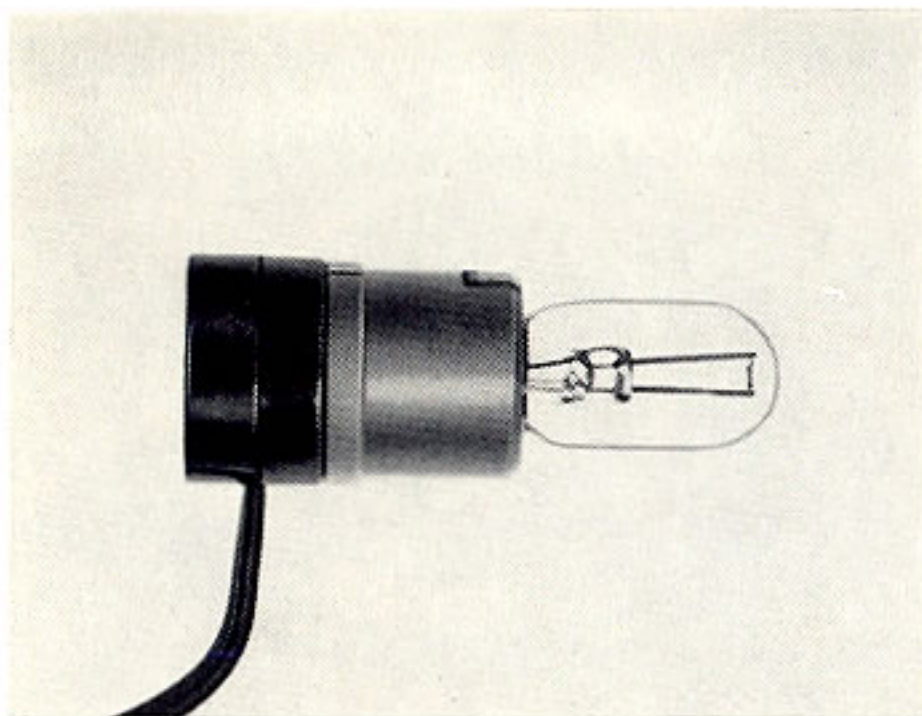
Changing bulb

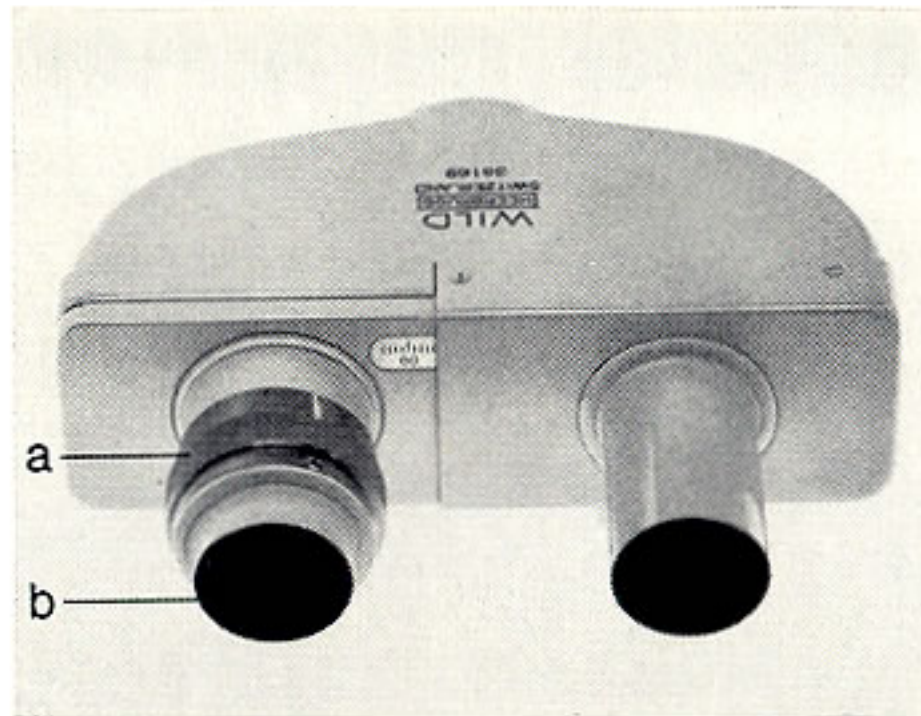Withdraw lamp socket from Köhler illuminator. Press bulb against socket and remove from bayonet by slight turn to left. Fit new bulb, replace lamp socket and secure clamping ring.

Adjusting binocular tube

Inter-pupillary distance: Slide left eyepiece (a) sideways until a single circular field is seen.

Anisometropia: Looking through right eyepiece with right eye, focus on a fine structure within specimen. Then look through left eyepiece with left eye and bring same structure into focus by turning diopter ring (b), without using fine adjustment a second time.

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Care of the microscope

Protect the instrument from dust, dampness, and acid or alkali vapours.
Do not store chemicals near the instrument.
Replace the microscope in its cabinet after use.
Use the dust cover if the microscope is left on the bench.
Clean the stand and the lens surfaces with a soft dusting brush from time to time.
Never use force when screwing the objectives in and out of the nosepiece.
Never touch the lenses, mirror or bulbs with the fingers.
Clean dirty lenses with a soft rag, moistened with Xylol.
Prevent dust from entering the tube (leave the eyepieces in the tube).
Never, on any account, unscrew the objectives, eyepieces or mechanical components.
Do not grease or oil the mechanical parts.
If the instrument no longer functions satisfactorily ask the advice of a specialist or your local WILD representative.

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Packing

If the instrument has to be re-packed for transportation, the objectives are unscrewed from the nosepiece, the eyepieces removed from the tubes, and all stored in the optics holder. The mirror or socket lamp can remain in the microscope base.

Important: The optics holder is fitted between the base and the stage, in order to protect the fine focussing mechanism. The stage must be manually lifted slightly in order to store the holder (see picture on page 3). Lower microscope body completely by means of coarse focussing control (1). Place metal hood over microscope and secure it using the two clips.

Care! If the microscope has a mechanical stage, its co-ordinates should be brought into an approximately central position.

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Wild Heerbrugg Limited